Reasons to choose the MASTER

1 Water Resistant



The concept of a washable handheld analog unit is measuring, the instrument can be cleaned with running water under a faucet The new structuraldesign of these units allows them to be (dust tight and water resistant).

washed whenever necessary. Rated IP65



2 Hygienic Design



The MASTER series was also designed with a smooth and easy-to-clean grip, eliminating the possibility of food and samples being trapped which can result in contamination and bacterial

3 Visibility



Efficiently using an analog unit depends on how easily the measured value can be read. Extensive research and the use of quality optics result in a measurement scale that is very bright, creating a high contrast to see clearly where the boundary intersects the memory lines

4 Automatic Temperature Compensation (ATC) 0.T.H.535



The newly designed ATC system allows the MASTER series to read more accurately than any comparable unit. ATAGO has applied for a patent on the design of the improved ATC system. The MASTER series reads more accurately than any comparable unit in the

CV-H-535-500 5 Durability



The MASTER series has passed all water resistance, dust resistance, and drop tests. instruments were subjected to water jets form four directions, dropped from a height of one meter onto an oak platfoam, and withstood the pressure changes of air travel.

6 Calibration



The α and H types can now be calibrated with the new waterproof calibration screw. All MASTER models can be calibrated by end-

7 Automatic Sample Distribution (ASD)



ASD ATAGO has developed a spoon-shaped tip that automatically spreads the sample across the prism. This makes the instrument even more user-friendly. saving valuable time.

8 Repair and Service



ATAGO has established a close network of customers by providing excellent customer service and support . When an ATAGO instrument needs servicing, not only is each customer treated with respect, but they are also provided with detailed methods and userspecific suggestions on sampling and taking measurements, based on individual

Reputation for Quality



ATAGO has built a firm reputation for reliability and quality during its over 70 years history. We are a strong company that exports to over 154 countries with continued growth expectations. ISO 9001 certified, along with our integrated system for producing and marketing our products, based on the "Plan-Do-Check-Act" cycle, is essential to the continued quality and cycle, is essential to the continued quality and to the promoting of the ATAGO name.


Simple Operation



1 Apply 1-2 drops onto the prism.



2 Close the daylight plate. View the scale through the eyepiece.

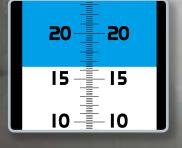

3 Read the measurement value where the boundary line intersects the scale.

Applications

-Measuring sugar concentration, concentration, Brix and Refractive Index-

Honey



Jam, Syrup



Jam, syrup, confectioneries

Soymilk



Condiments



Condiment, seasonings, ketchup, various sauces, sov sauce

Wine



Grapes, wine must

Fruit juice, soft drink, carbonated beverage

Drink

▶ Cooking



Chinese noodle soup, various soup, various sauces

Agriculture



Fruit and vegetable juice, leaf juice, stem juice, root juice

Salinity



sea water, pickles

Machine Tools



Firefighting foam, soluble cutting oil

Chemicals and medicine



Brine, deicing solution, medicine, chemical, cosmetics, cleaning solution

Clinical



Urine, serum

Optional Accessories



Daylight Plate Parts.No. RE-2315-60M Parts.No. RE-2315-61M

Small Volume Daylight Plate With Hinge Pin







Sucrose solution (for Brix confirmation)

Parts.No. RE-110010 10% Sucrose Parts.No. RE-110020 20% Sucrose Parts.No. RE-110030 30% Sucrose Parts.No. RE-110040 40% Sucrose Parts, No. RE-110050 50% Sucrose Parts.No. RE-110060 60% Sucrose

All ATAGO refractometers are designed and manufactured in Japan.



Headquarters: The Front Tower Shiba Koen, 23rd Floor 2-6-3 Shiba-koen, Minato-ku, Tokyo 105-0011, Japan TEL: 81-3-3431-1943 FAX: 81-3-3431-1945







ATAGO RUSSIA Ltd.

ATAGO BRASIL Ltda.

ATAGO ITALIA s.r.l.

ATAGO NIGERIA Scientific Co., Ltd.

ATAGO CHINA Guangzhou Co., Ltd.

TEL: 66-21948727-9 TEL: 55 16 3913-8400 TEL: 39 02 36557267

TEL: 1-425-637-2107


TEL: 86-20-38108256 TEL: 7-812-777-96-96

TEL: 91-22-28544915 40713232 customerservice@atago-india.com customerservice@atago-thailand.com customerservice@atago-italia.com

> info@atago-china.com info@atago-russia.com

* Specifications and appearance are subject to change without notice.

Copyright © 2015 ATAGO CO., LTD. All rights reserved. ENV.06 15091000PP Printed in Japan

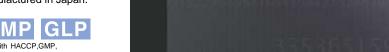













The Ultimate Hand-Held Refractometer Unsurpassed

in Function, Operation, and Design

MASTER





MASTER series

Choose the right model for your sample, based on the features and materials of the instrument.



Features 6 types

Both the α and T types are equipped with the Automatic Temperature Compensation feature. Choose the α type for the added feature of IP 65 water-resistance. The H type is the heat-resistant version of the α type.

Series / Model	а	Т	M	Н	53S	500
Water resistant		_	_	•		
Automatic temperature compensation	•	•	_	•	•	_
Heat resistant	_	_	_	•	- \	•
Milky sample	_	_	_	_	•	_

MASTER-53S

for Milky sample

0-90% Wide Brix Range

Opaque, inhomogeneous liquids can be difficult to get a clear boundary line with a conventional refractometer. This model is recommended for cream-based sauces, yogurt, mayonnaise, and various emulsions used in industrial processing.








■ Understanding model names

A model name consists of the scale (sample name), material, and feature type.



Material 2 types Choose either metal or plastic.

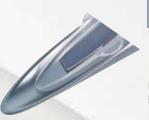

Die-cast aluminum

Ideal for the agriculture industry, metal working fluids, and various high-temperature samples. Resistant to organic solvents.



Plastic

PBT (polybutylene terephthalate)

Resistant to salty and acidic samples.

■ Brix of Various Samples Refractive index and Brix of various types of samples

Brix	(%)	Food industry			Chemical industry	Medical Science industry			_	(Brix ory No.	
(Ri)	^	Drinks				Methanol					
1.333	0	Green tea Beer	Fruits /	Seasonings Miso soup	Cutting oil (emulsion)	Saline Aloe extract Sea water	(1) 0 5	10 0 5	20 0 5 20 %	53 0 5	500 0 5
1.34 —	5	Coffee	Tomato Strawberry Lemon		Nickel plate liquid	Chlorella extract	33 %	10 %	20 %	53 %	90 %
1.35 —	10	Soy milk Cola Milk	Orange Apple Prince melon	Beef curry	Gelatin liquid Freon113	Serum					
	15 ·			Demiglace (sauce) Orange jelly							
1.36 -	20	Lactic acid- bacteria be		Canned syrup	Silicone solution (emulsion)	Acetone Ethanol					
1.37 —	30			Gailliod Sylap	Electrolyte Freon 11	Acetic acid Saturated saline		₿			
1.39 —	30			Ketchup Soy sauce Jam	Acrylonitrile			2 28 5			ı
1.40 -	40			Sauce A Low-sugar jam				62			
	50 ··			Egg yolk	Silicone oil	Octane		70	. 4		
	60	Fats Butter			Ethylene glycol Propylene glycol		3		45 \$ 82		
	70	Olive oil		Honey Jam	Diesel oil		58 \$		%		
1.50	80			Honey Starch Syrup	Ricinus	Glycerin	90				
1.50	90				Polyvinyl chloride (PCV) Styrene Asphalt						

■ Specifications

Brix Scale

Range	Series	(z c	1	Γ	Ņ	Л
(Minimum)	Material	Metal	Plastic	Metal	Plastic	Metal	Plastic
4	Model	MASTER-α	MASTER-Pα	MASTER-T	MASTER-PT	MASTER-M	MASTER-PM
0.000/	Cat.No.	2311	2391	2312	2392	2313	2393
0 to 33%	Accuracy	±0.2% *±0.1	(10 to 30°C)	±0.2% *±0.1	% (10 to 30°C)	±0.2%	*±0.1%
(0.2%)	Size & Weight	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g

2	Model	MASTER-2α	MASTER-2Pα	MASTER-2T	MASTER-2PT	MASTER-2M	MASTER-2PM
201.000/	Cat.No.	2321	2941	2322	2942	2323	2943
28 to 62%	Accuracy	±0.2% *±0.1% (10 to 40°C)		±0.2% *±0.1% (10 to 40°C)		±0.2% *±0.1%	
(0.2%)	Size & Weight	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g

2	Model	MASTER-3α	MASTER-3Pα	MASTER-3T	MASTER-3PT	MASTER-3M	MASTER-3PM
501.000/	Cat.No.	2331	2951	2332	2952	2333	2953
58 to 90%	Accuracy	±0.2% *±0.1% (10 to 40°C)		±0.2% *±0.1% (10 to 40°C)		±0.2%	*±0.1%
(0.2%)	Size & Weight	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90q	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90q

4	Model	MASTER-4 a	MASTER-4Pα	MASTER-4T	MASTER-4PT	MASTER-4M	MASTER-4PM
45. 000/	Cat.No.	2341	2961	2342	2962	2343	2963
45 to 82%	Accuracy	±0.2% *±0.1% (10 to 40°C)		±0.2% *±0.1% (10 to 40°C)		±0.2% *±0.1%	
(0.2%)	Size & Weight	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g

40	Model	MASTER-10α	MASTER-10P α	MASTER-10T	MASTER-10PT	MASTER-10M	MASTER-10PM
10	Cat.No.	2371	2981	2372	2982	2373	2983
0 to 10%	Accuracy	±0.2% *±0.1	% (10 to 30°C)	±0.2% *±0.1	% (10 to 30°C)	±0.2%	*±0.1%
(0.1%)	Size & Weight	3.2×3.4×20.3cm, 155g	3.2×3.4×20.3cm, 105g	3.2×3.4×20.3cm, 155g	3.2×3.4×20.3cm, 105q	3.2×3.4×20.3cm, 155q	3.2×3.4×20.3cm, 105q

-	20	Model	MASTER-20 α	MASTER-20P α	MASTER-20T	MASTER-20PT	MASTER-20M	MASTER-20PM
	20	Cat.No.	2381	2991	2382	2992	2383	2993
	0 to 20%	Accuracy	±0.2% *±0.1% (10 to 30°C)		±0.2% *±0.1% (10 to 30°C)		±0.2%	*±0.1%
- 1	(0.1%)	Size & Weight	3.2×3.4×20.7cm, 165g	3.2×3.4×20.7cm, 110g	3.2×3.4×20.7cm, 165g	3.2×3.4×20.7cm, 110g	3.2×3.4×20.7cm, 165g	3.2×3.4×20.7cm, 110g

52	Model	MASTER-53α	MASTER-53Pα	MASTER-53T	MASTER-53PT	MASTER-53M	MASTER-53PM
53	Cat.No.	2351	2971	2352	2972	2353	2973
0 to 53%	Accuracy	±0.5% *±0.25	% (10 to 30°C)	±0.5% *±0.25	% (10 to 30°C)	±0.5%	*±0.25%
(0.5%)	Size & Weight	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g	3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g

► Heat resistant

Series		Н							
Material			Metal						
Model	MASTER-50H	MASTER-80H	MASTER-93H	MASTER-100H					
Cat.No.	2354	2364	2374	2384					
Range	Brix 0.0 to 50.0%	Brix 30.0 to 80.0%	Brix 45.0 to 93.0%	Brix 60.0 to 100.0%					
Minimum			0.5%						
Accuracy		±0.5% *±0.25%							
Size & Weight		3.2×3.4	× 16.8cm, 130g						

► Milky sample ► Wide Brix

Series	53S	500		
Material	Me	etal		
Model	MASTER-53S	MASTER-500		
Cat.No.	2355	2363		
Range	Brix 0.0 to 53.0%	Brix 0.0 to 90.0%		
Minimum	0.5%	1.0%		
Accuracy	±0.5% *±0.25%	±1% *±0.5%		
Size & Weight	3.2×3.4× 16	.8cm, 130g		

Special Scale

Sample	Coolant	Battery Coolant	Battery Coolant	Refractive index	Honey
Series	M	M	M	M	α
Material	Plastic	Plastic	Plastic	Metal	Plastic
Model	MASTER-BR	MASTER-BC	MASTER-BCF	MASTER-RI	MASTER-HONEY
Cat.No.	2930	2931	2932	2612	2524
Scale ① propylene glycol ② ethylene glycol		ethylene glycol freezing temp propylene glycol freezing temp sterp duid d20/20	ethylene glycol freezing temp propylene glycol freezing temp stery fluid d20/20	Refractive index	Honey moisture
Range	① 0 to 70% 0 to -50℃ ② 0 to 70% 0 to -50℃	① 0 to -50℃ ② 0 to -50℃ ③ 1.150 to 1.300	① 32 to -60°F ② 32 to -50°F ③ 1.150 to 1.300	1.435 to 1.520	12.0 to 30.0%
finimum Scale	5% 5℃	① 5℃ ② 5℃ ③ 0.01	① 10°F ② 10°F ③ 0.01	0.001	0.2%
Size & Weight		3.2×3.4×16.8cm, 90g		3.2×3.4×16.8cm, 130g	3.2×3.4×16.8cm, 90g

Sample	Seawater	Seawater	Salinity	Salinity	Salinity (Wide range)	Salinity (Wide range)
Series	α	M	α	M	α	M
Material	Plastic	Plastic	Plastic	Plastic	Plastic	Plastic
Model	MASTER-S/Mill α	MASTER-S/Mill M	MASTER-S10 α	MASTER-S10 M	MASTER-S28α	MASTER-S28 M
Cat.No.	2491	2493	2471	2473	2481	2483
Scale	Salinity ② Specific gravity		Sodium chloride		Sodium chloride	
Range	① 0 to 100‰ ② 1.000 to 1.070		0.0 to 10.0g/100g		0.0 to 28.0g/100g	
linimum Scale	① 1‰ ② 0.001		0.1g/100g		0.2g/100g	
Measurement Accuracy	① ±2‰ *±1‰ ② ±0.001 *±0.0005 (10 to 30°C)	_	±0.2g/100g *±0.1g/100g (10 to 30°C)	_	±0.2g/100g *±0.1g/100g (10 to 30°C)	_
International rotection Class	IP65		IP65	_	IP65	_
Size & Weight	3.2×3.4×20.7cm, 110g		3.2×3.4×20.3cm, 105g		3.2×3.4×16.8cm, 90g	

IP65 = dust-tight and protected against water jets (except eye piece)