



Powerful Performance Comes in Small Packages





### ATAGO's new polarimeter offers robust performance and ultimate efficiency in small footprint.

Thanks to improvements made in structural design, the POL-1/2 needs only half the space of a conventional model. The space-saving design promotes installation flexibility. A lightning-fast response time of 60° per second optimizes operational efficiency.

#### Looks can be deceiving.

Though it is half the size of a traditional polarimeter, the POL- ½ provides better resolution than most.

There is no time to waste.

The response time is 60° per second. That is 15 times faster than before.

It cannot be easier. The LCD touchscreen color display makes all operations fast and easy. Data transfer is simple with the use of a USB flash drive. The sample temperature is measured directly and accurately.

Always aim high.

With new optical technology, the world's highest class resolution of 1/10000° has been achieved.

Get validated. Stay compliant.

We provide Installation Operational Performance Qualification (IQ/OQ) documentation and support to ensure compliance with all appropriate regulations.

## **Powerful Performance Comes in Small Packages**



POL - 1/2

POL-1/2 Cat.No.5271

Pharmaceutical

Various medications and vitamins

## Concept

## With a 100mm cell, measure at an accuracy of $\pm 0.002^{\circ*}$ and resolution of 0.0001°.

\* Less than 1° angle of rotation

Compliance with pharmacopeias and other regulatory requirements worldwide is within easy grasp.





## **Temperature Control**

### The optional Peltier temperature control unit delivers perfectly temperature-controlled measurements.


This Peltier temperature control unit does not require the use of water. Enjoy stress-free temperature control without worrying about a water leak or cleaning.



POL-1/2 with Peltier temperature control Cat.No.5272



Target temperature

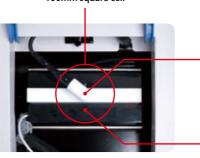


100mm square cell

Temperature sensor for sample (standard accessory) Choose either the temperature of the cell holder or the temperature of the sample liquid itself to be measured.

Advantage of using a square cell The thermostat is in contact with the bottom and both sides of the cell, enabling quick and accurate temperature control.

#### Measurement results when using a Peltier temperature control unit - The closer the sample temperature is to the target temperature, the more quickly the stability is obtained.


Specific rotation of saccharose Angle of rotation of water cell: 100mr 67.15 21.05 0.0015 21.03 - 67.10 0.0010 67.05 21.00 + [α] 0 0005 h M M M 67.00 20.98 0.0000 66.95 20.95 -0.0005 66.90 20.93 -0.0010 -0.0015 66.85 20.90 50 50 10 40 10 20 30 40 60 20 30 60 0 Time (mir Time (min) Water Saccharose (surcorse)

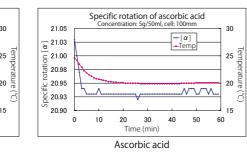

The hollow-shaft motor turns the analyzer directly. Measure at 60° per second, the best response speed in its class.

#### LED light source \*Only 589nm is available


explicitly listed in traditional pharmacopeias, it is regarded as compliant in a practical sense for the with the results obtained with other measuring devices listed in pharmacopeias.

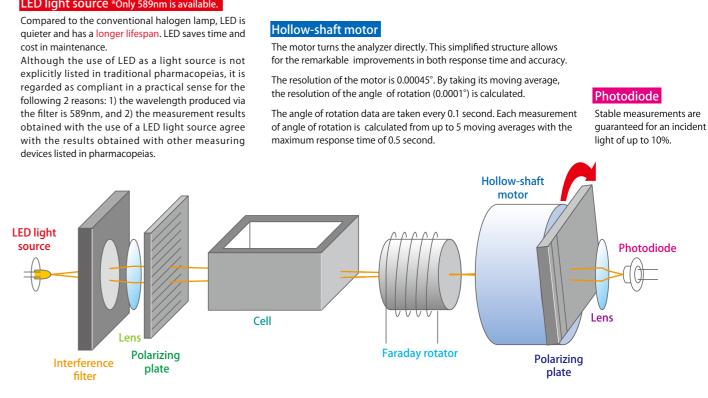

Polarizing plate A polarizer with a high extinction ratio is used

#### Interference filter

A high performance filter with a FWHM of 2.5nm is used.

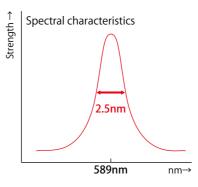


# Cell

## **Mechanism**

### The hollow-shaft motor delivers industry's fastest measurement.

In addition to the standard square shape, cylindrical cells are also available upon request. Choose from 10mm, 20mm, 50mm, or 100mm. The exact cell length is entered in the unit

for measurement. \*The exact optical path length cannot be entered in the unit. The volume capacity of a standard cell with an optical path length of 100mm is approximately 12ml. The smallest cell is 10mm in optical path length and 0.1ml in volume. Cells are available in various volume capacities of 0.1 to 30ml to meet many different application needs.

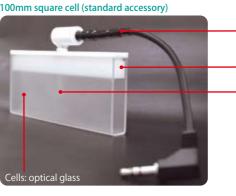

Temperature sensor for sample (standard accessory)

Lid: PTFE


Temperature sensor: PFA The sensor probe is directly submerged in the sample liquid for optimal accuracy.

## Function

### The color LCD touchscreen delivers quick and smooth operations.


With the intuitive user interface, it is extremely easy to view measurement results and adjust settings.


#### Measurement result screens

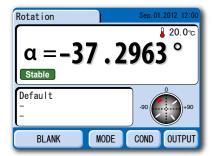


#### Settings and data input screens

|--|

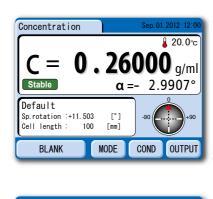

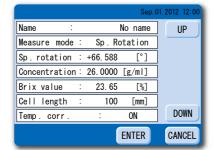

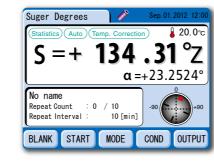

		Sep.01.2012 12:00
1 ~ 5 6 ~ 10 1 :Default 2 : 3 : 4 : 5 :	Name : Mode : Sp. rotation Concentration Brix Cell length Temp corr . Corr . coef . Statistics Auto measure Repeat number Auto interval Responce Wave length	: 23.65 [%] : 100 [mm] : 0FF
	S	AVE CANCEL


#### Example of continuous measurement settings (mutarotation of glucose)

Examples of setting items Auto measure · ON Repeat number: 99 Auto interval: 4 Response: 5 (default)

Data storage and export

[°] [°C]

Mutarotation of a substance can be measured by setting the number of measurements and time interval between measurements in the continuous measurement mode

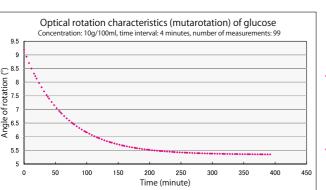

#### Data storage

The graph on the left was created in Excel from the text file of the measurement results transferred to a PC via a USB flash drive.

#### Protection of settings

Up to 10 sets of settings can be stored and protected with passwords.

> ATAGO POL-1/2 2012 09/01 12:00 NAME : Test\_Sample\_01 LAMP : 589 [rm] CONCENTRATION : 9,26000[g/m] o(n−1)= 0.0012 (\*) C.V. = 0.0031 [½]

> > (up to 10 measure

Cells



#### DP-AD24 dot matrix printer



Prints out measurement data. Data can be exported to either this printer or a USB flash drive.



\* Connection of circulating constant temperature water bath The circulation flow should be set to 12 L / min as a guide, and the maximum head pressure should not exceed 5 m. Prepare rubber piping with 11 mm O.D. / 7 mm I.D.

#### Product Name Part No. RE-86010 Square type cell hol RE-82100 Square cell 100 RE-82101 Square cell 50 RE-82102 Square cell 20 RE-82103 Square cell 10 Square cell lid 100 RE-88100 RF-88101 Square cell lid 50 RE-88102 Square cell lid 20 Square cell lid 10 RE-88103 Plug for cell lid, 2pc RE-88200 RE-89100 Temperature sensor RE-87000 Cell holding parts RE-82104 Square micro cell 10 Square micro cell 50 RE-82105 RF-86103 Aperture RE-86011 V-type cell holder Cylindrical cell 100 RE-88014 RE-88015 Cylindrical cell 50 RE-88016

#### Cylindrical cell 20 RE-88017 Cylindrical cell 10 RE-88108 Round cell lid RE-82106 Jacket cell A1 RE-82107 Jacket cell A3 RE-82108 Jacket cell B1 RE-82109 Jacket cell C1 RE-82110 Jacket cell C2 RE-82111 Jacket cell C3 Jacket cell C4 RE-82112 Cat.No.3133 DP-AD24 RE-89403 Standard roll paper

Ink ribbon for standa

TCS-1

Angle of rotation

ROTATION « = +10.0193 - 16.4

ATAGO POL-1/2 2012 09/01 12:00 NAME : Test\_Sample\_01 LAMP : 589 [nm] CELL LENGTH : 100 [mm]

Specific rotation

ROTATION

ATAGO POL-1/2 2012 03/01 12:00 NAME : Test\_Sample\_01 LAMP : 589 [nm] CONCENTRATION : 0.26000[g/ml] CELL LENGTH : 100 [mm]

SPECIFIC ROTATION [4] = +38.533 [°]

= +10.0186 [°] TEMP = 16.4 [°C]

Angle of rotation, specific rotation, concentration, International Sugar Scale readings as well as

statistical computation results can be exported. Measurement results can be either saved as a

text file to a USB flash drive or printed to a printer (sold separately).

USB flash drive (1GB/FAT16) is standard accessory.

SPECIFIC ROTATION (d) = +38.484 [°] ROTATION d +10.0191 16.4 [°] [°C] TEMP = Specific rotation (temperature compensated)

ATASO POL-1/2 2012 09/01 12:00 NAME : Test\_Sample\_01 LAMP : SS9 [nm] CONCENTRATION : 0.26000[g/ml] CELL LENGTH : 100 [nm] TEMP CORR. COEF. : 0.00037

	SPECIFIC COUNT 1 2 3 4 5 6	DF +38.5 +38.5 +38.5 +38.5	ATA 532 530 531 532 530	C() (*) TEMP 16.3 16.3 16.3 16.3 16.3 16.3
	MEAN = σ(n-1)= C.V. =	+38. 0.0	531 0010	[°] [°] [%]
-	Average &	stanc	lard d	leviatior

ATAGO POL-1/2 2012 03/01 12:00 NAME : Test.Sample.01 LAMP : S80 [rmn] CONCENTRATION : 0.26000(g/ml] CELL LENGTH : 100 [mm1]

SPECIFI	C ROTATION	
COUNT	DATA	
1	+38.530	16.3
2	+38,538 +38,528 +38,528 +38,528 +38,528 +38,528 +38,528 +38,528 +38,528 +38,528 +38,528 +38,528	16.3
3	+38.528	16.3
4	+38.528	16.3
5	+38.528	16.3
6	+38.526	16.3
7	+38.530	16.3
8	+38.528	16.3
9	+38.528	16.3
10	+38.527	16.3

Statistical computation results

RE-89402 RE-86012 Cat.No.1922 60-C4

## Accessories

## A wide variety of cell options, printer deliver optimized measurement and data processing solutions.


	Detail
lder	For square cell, without cell lid, applicable for circulated water at constant temperature (same as standard supplied holder)
	Optical path length 100 mm, sample amount 12 mL, without cell lid (spare)
	Optical path length 50 mm, sample amount 6 mL, without cell lid
	Optical path length 20 mm, sample amount 2.4 mL, without cell lid
	Optical path length 10 mm, sample amount 1.2 mL, without cell lid
	For square cell 100, made of fluorine resin, with a hole for temp sensor probe
	For square cell 50, made of fluorine resin, with a hole for temp sensor probe
	For square cell 20, made of fluorine resin, with a hole for temp sensor probe
	For square cell 10, made of fluorine resin, with a hole for temp sensor probe
CS .	For cylindrical and square cells, spare
r for sample	For sample temperature measurement, standard accessory (spare)
	For sample temperature measurement, using 10 mm / 20 mm square cell
00	Optical path length 100mm, sample amount 2mL, cell lid included, the aperture for square micro cell required
0	Optical path length 50mm, sample amount 1mL, cell lid included, the aperture for square micro cell required
	For square micro cell
	For cylindrical and jacket cells
	Optical path length 100mm, sample amount 29.8mL, without cell lid, V-type cell holder required
	Optical path length 50mm, sample amount 14.9mL, without cell lid, V-type cell holder required
	Optical path length 20mm, sample amount 6.0mL, without cell lid, V-type cell holder required
	Optical path length 10mm, sample amount 3.0mL, without cell lid, V-type cell holder required
	For cylindrical, jacket, and square micro cells, made of fluorine resin, with a hole for temperature sensor probe
	Optical path length 100mm, sample amount 6.4mL, V-type cell holder required
	Optical path length 20mm, sample amount 1.3mL, V-type cell holder required
	Optical path length 100mm, sample amount 2mL, V-type cell holder required
	Optical path length 100mm, sample amount 1mL, V-type cell holder required
	Optical path length 50mm, sample amount 0.5mL, V-type cell holder required
	Optical path length 20mm, sample amount 0.2mL, V-type cell holder required
	Optical path length 10mm, sample amount 0.1mL, V-type cell holder required
	Plain paper printer - Connection cable included (0.75m)
	4 rolls in a set
ard paper printer	1 piece
	Peltier temperature control unit
	Circulating constant temperature bath (4L / min)

# Specifications

#### Polarimeter

#### Dimensions

0000000000 250  $2 \times \phi 8.2 <$ 

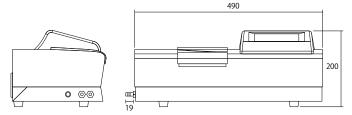

#### Technical specifications

Light source Polarizer Wavelength Detector Observation tube length Measurement range Accuracy


Repeatability Resolution Response time Measuring principle Thermometer location Temperature range Display Input method Measurement item

Data processing Data storage Network connection Power supply Power consumption Dimensions Weight

LED Polarizing film 589nm Photodiode Up to 100mm  $\pm$  90° angle of rotation  $\pm 0.002^{\circ}$  of displayed value (less than 1°), relative precision of  $\pm 0.2\%$  (greater than or equal to 1°)  $+0.002^{\circ}$ 0.0001° 60°/second max. Symmetric angle oscillation optical null method At the cell holder or inside the cell 0.0 to 99.9℃ 5.7-inch color LCD Touchscreen Angle of rotation, specific rotation, International Sugar Scale, concentration, purity Continuous measurement, statistical computation Printer (sold separately), USB flash drive Ethernet AC100V to AC240V, (50/60Hz) 100VA 490 (W) x 250 (D) x 200 (H) mm 13kg

### Peltier temperature control unit

#### Dimensions

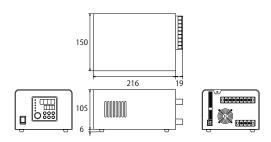

#### Technical specifications

Temperature control method	Auto-tuning PID temperature controller
Temperature range	15℃ to 35℃
Control accuracy	$\pm$ 0.1°C (resistance thermometer)
Display accuracy	0.1℃
Resolution	0.1℃
Ambient temperature range	0 to 40℃
Ambient humidity range	Less than 90% RH, no condensation
Safety features	Fuse (interrupts excessive current), buzzer alarm
Power supply	AC100V to AC240V, (50/60Hz)
Power consumption	150VA
Dimensions*	150 (W) x 216 (D) x 105 (H) mm
Weight	Approx. 2.5kg

\* Rubber feet, terminal blocks, and any other protuberances are excluded.

All ATAGO polarimeters are designed and manufactured in Japan.

### ATAGO CO., LTD.

Headquarters: The Front Tower Shiba Koen, 23rd Floor C-6-3 Shiba-Koen, Minato-ku, Tokyo 105-0011, Japan TEL : 81-3-3431-1943 FAX : 81-3-3431-1945 overseas@atago.net http ://www.atago.net/ TEL: 1-425-637-2107 customerservice@atago-usa.com

TAGO INDIA hetruments Pvt. Ltd. TEL : 91-22-2833-8038 / 8076 customerservice@atago-india.com

CATAGO (THAILAND) Co., Ltd. TEL : 662-982-8718-9 customerservice@atago-thailand.com

TEL : 55 16 3916-6000 customerservice@atago-brasil.com

ATAGO ITALIA s.r.l. TEL : 39 2 36557267 customerservice@atago-italia.com

ATAGO CHINA Guangzhou Co., Ltd. TEL : 86 20-38106065 info@atago-china.com





\*Specifications and appearance are subject to change without notice.