

Sugar-Free Drinks

Diet Drinks

etc.

Specially developed for very low concentration solutions, Brix 8% or less!!

Hand-held Refractometer
(For very low concentration samples)

N-8α

Cat.No.2360



Applications

Food industry :

- Sugar-Free
- Diet drink
- very thin soup

etc.

Measuring Method




①Place a couple of drops of sample onto the prism.



②Close the day-light plate gently.



③Look through the eyepiece and measure the value.

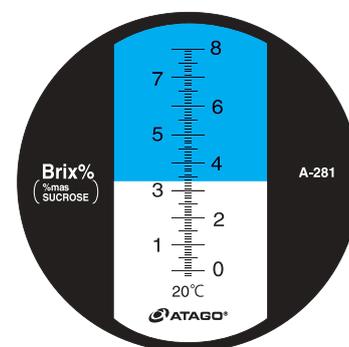

Features

- Bright, High Contrast scale view
- Scale memory lines are wider for cleaner reading of the boundary line
- Requires just one or two drops of sample
- Quick & easy measurements

Specifications

Measurement range: Brix0.0 to 8.0%
Minimum scale : 0.1%
Dimensions : 4(W)×4(D)×21(H)cm
Weight : 260g

All ATAGO refractometers are designed and manufactured in Japan.

ATAGO CO.,LTD.

<http://www.atago.net/> overseas@atago.net

Headquarters: The Front Tower Shiba Koen, 23rd Floor
2-6-3 Shiba-koen, Minato-ku, Tokyo 105-0011, Japan
TEL : 81-3-3431-1943 FAX : 81-3-3431-1945



HACCP GMP GLP

ATAGO products comply with HACCP,GMP, and GLP system standards.

- **ATAGO U.S.A., Inc.**
- **ATAGO INDIA Instruments Pvt.Ltd.**
- **ATAGO THAILAND Co.,Ltd.**
- **ATAGO BRASIL Ltda.**
- **ATAGO ITALIA s.r.l.**
- **ATAGO CHINA Guangzhou Co.,Ltd.**
- **ATAGO RUSSIA Ltd.**
- **ATAGO NIGERIA Scientific Co.,Ltd.**

- TEL : 1-425-637-2107 customerservice@atago-usa.com
- TEL : 91-22-28544915, 40713232 customerservice@atago-india.com
- TEL : 66-21948727-9 customerservice@atago-thailand.com
- TEL : 55 16 3913-8400 customerservice@atago-brasil.com
- TEL : 39 02 36557267 customerservice@atago-italia.com
- TEL : 86-20-38108256 info@atago-china.com
- TEL : 7-812-777-96-96 info@atago-russia.com
- TEL : 234-707-558-1552 atagonigeria@atago.net

* Specifications and appearance are subject to change without notice.