### **Optional Accessories**

Part No.	Part name		
< Container	>		
RE-79100	15mL Beaker		
RE-79101	100mL Beaker		
RE-78141	Cup Adapter (with 100pcs cups) * 50pcs of paper cups and 50pcs of plastic cups are included.		
RE-79102	Paper Cup (90mL, 100pcs)		
RE-79103	Plastic Cup (90mL, 100pcs)		
< Ultra Low	Adapter (ULA)>		
RE-77120	Ultra Low Adapter (ULA)-		
	Sample Adapter for Low Viscosity Sample		
	<ul> <li>Sample cylinder</li> </ul>	Cylinder holder	
	• Hook	Hook holder	
	<ul> <li>UL spindle</li> </ul>	• UL stand	
	Extension (threaded tip)		
RE-77107	UL spindle (with hook and hook holder)		
RE-77121	Sample cylinder (with cap and o-ring)		
RE-77117	UL spindle 3pcs (with hook and hook holder)		
< Temperati	ure sensor >		

Part No.	Part name	
< Spindle >		
RE-77104	A1 Spindle	
RE-77105	A2 Spindle	
RE-77106	A3 Spindle	
RE-77114	A1 Spindle 5pcs	
RE-77115	A2 Spindle 5pcs	
RE-77116	A3 Spindle 5pcs	
RE-77100	Set of spindles (A1,A2,A3)	
< Viscosity S	Standard Liquid >	
RE-89030	Viscosity Standard Liquid 2	100mL
RE-89031	Viscosity Standard Liquid 5	100mL
RE-89036	Viscosity Standard Liquid 200	100mL
RE-89037	Viscosity Standard Liquid 500	100mL
RE-89038	Viscosity Standard Liquid 1000	100mL
RE-89039	Viscosity Standard Liquid 2000	100mL

\* Standard liquid with JCSS calibration certificates are available (JS2.5 to JS160000). Contact ATAGO for further details.

RE-75540 Temperature sensor

#### Specifications Measurement range 1 to 350,000,000mPa·s, 1 to 350,000,000cP

Model	VISCOTM	VISCO™-895
Cat.No.	6800	6820
Materials	Housing: SUS, Aluminum · Legs, and stand + screw: SUS	Housing, legs, and stand + screw: Aluminum
Dimensions and Weight	12×12×20cm, 1.2kg (main unit only), Stand+screw : 0.5kg	12×12×20cm, 895g (main unit only), Stand+screw : 275g
	Small volume beaker attachment:	Small volume beaker attachment:
	0.1kg	0.1kg

#### The body, legs and stage of the VISCO<sup>™</sup>-895 are made from light-weight aluminum.



Contents	
• Main unit	1
· Stand:	1
· S Beaker (15mL)	1
· L Beaker (100mL)	1
· AC adapter	1
· Spindles (A1, A2 and A3)	one each
· Temperature sensor	1
· Small volume beaker attachmer	nt 1
· USB Mini-B cable (1m)	1
· 1.5V AA alkaline batteries	4
<ul> <li>Instruction manual</li> </ul>	1
<ul> <li>Inspection certificate</li> </ul>	1
· Spindle stand	1
· Protective cap	1
· Carrying case	1

TEL: 1-425-637-2107

Measurement	Viscosity · Temperature · Torque%			
Scales	viscosity · temperature · torque/0			
Measurement	Viscosity	A1 50 to 200,000mPa·s, 50 to 200,000cP		
Range		A2 100 to 600,000mPa·s, 100 to 600,000cP		
		A3 500 to 2,000,000mPa·s, 500 to 2,000,000cP		
		(1mPa·s=1cP)		
	Torque	0.0 to 100.0%		
		(recommended torque : 10.0 to 100.0%)		
	Temperature	0.0 to 100.0°C / 32.0 to 212.0°F		
Resolution	Viscosity	lower than 100mPa·s : 0.01mPa·s		
		100mPa·s or higher lower than 10,000mPa·s : 0.1mPa·s		
		10,000mPa·s or higher : 1mPa·s		
	Torque	Lower than 10% : 0.01%		
		10% or higher : 0.1%		
	Temperature	0.1°C / 0.1°F		
Measurement	Viscosity	±1% (Full scale)		
Accuracy	Temperature	±0.2°C/±0.4°F		
Speed	0.5 to 250 rp	m, Number of speeds : 20		
Sample	10.0 to 40.0%	C / 50.0 to 104.0°F		
Temperature Range	10.0 to 40.0	07 00.0 10 104.0 1		
Ambient	10 to 40°C			
Temperature	10 10 40 0			
Computer Output	Output : USB - PC			
Battery Life (Approx.)	Approx. 7 ho	urs (continuous operation at 60rpm)		
Power Supply	DC6V (AA alk	aline batteries 1.5V×4)		
	AC adapter : AC100 to 240V, 50/60Hz			

All ATAGO products are designed and manufactured in Japan.



https://www.atago.net/ overseas@atago.net

ATAGO INDIA Instruments Pvt, Ltd, CATAGO BRASIL Ltda. CATAGO ITALIA s.r.I. CATAGO CHINA Guangzhou Co., Ltd. TEL : 86-20-38108256 info@atago-china.com CATAGO NIGERIA Scientific Co., Ltd. TEL : 234-707-558-1552

TEL: 91-22-28544915 / 40713232 customerservice@atago-india.com TEL: 66-21948727-9, 66-21171549 customerservice@atago-thailand.com TEL: 55 16 3913-8400 customerservice@atago-brasil.com TEL: 39 02 36557267 customerservice@atago-italia.com TEL: 7-812-777-96-96 info@atago-russia.com atagonigeria@atago.net CATAGO KAZAKHSTAN Ltd. TEL : 7-727-257-08-95 info@atago-kazakhstan.com

customerservice@atago-usa.com


\* Specifications and appearance are subject to change without notice.





Copyright © 2019 ATAGO CO., LTD. All rights reserved. ENV.07 19120300WA Printed in Japan



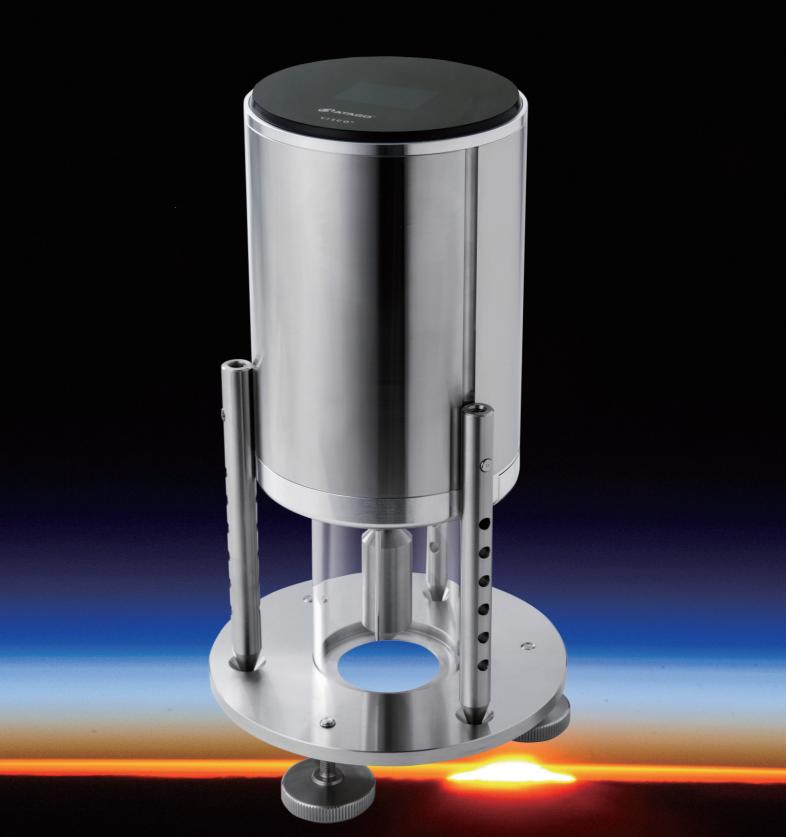







# ATAGO: Creating the Perfect Fusion of Innovation, Technology and Simplicity

## **ONE TOUCH™**, **ONE HAND™** and **ONE BUTTON™**.

Presenting the VISCO, a brand-new way of measuring viscosity with 3 simple **"ONE's."** 

# **ONE BUTTON**<sup>TM</sup>

Operation requires only one dial button. All operations can be performed with the simple act of "sliding" or "pushing" the dial button. No more accidental operations due to pushing the wrong button.

Measurement

# **ONE TOUCH™**


VISCO is very easy to set-up. The spindle can be attached with just "one touch" simply insert the spindle in the instrument. Absolutely no complicated set-up required.

Set-up

# **ONE HAND**<sup>™</sup>


Measurement preparation can easily be done with just one hand. Place the beaker underneath the pre-set area and place the instrument on the stand. No troublesome height adjustment necessary.

Preparation

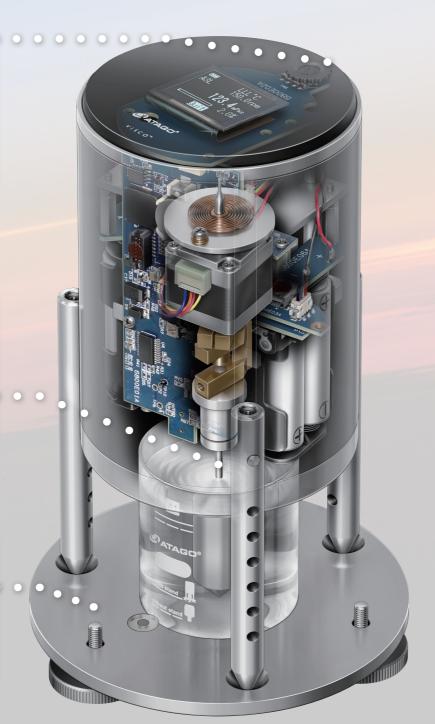


# Uses Only 1/33 of the Standard Sample Amount

Standard measurement instruments for viscosity require a large amount of sample (500mL). VISCO is capable of taking measurements with just 15mL of sample. This is roughly 1/33 of the standard sample amount. Measurement can be done with only a small amount of sample, resulting in less waste of valuable sample and a significant reduction in cost.

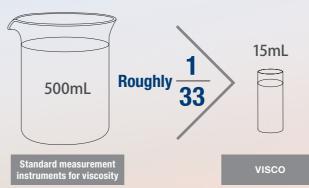

# Easy to Read, Fully Digital Display

A fully digital display allows for anyone to quickly and easily read results. The simple display is easily and readily understood.  $\therefore$  20.0°C  $\Rightarrow$  20.0°C  $\ddagger$ 

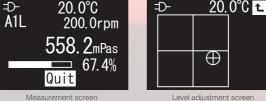

## Never Take Any Glassware (on site) Ever Again

Measurement can also be taken using paper cups other than those that come with the instrument. A disposable container requires no washing after each measurements. This makes it possible to safely measure even on site where no glassware is allowed.

• VISCO Package: Package A P.4


Disposable container can be purchased independently.



# **Compact and Easily Carried with One Hand**


VISCO's sleek dimensions and weight (main unit: 12x12x20cm, 1.2kg) make it compact and easily carried with one hand.

The instrument's legs can be folded up, making it further compact and allowing for even greater storage capabilities.

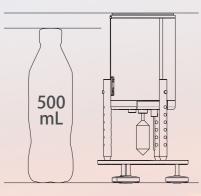

## **Quick Measurements Anywhere**

Not only does VISCO run on AC power, but it can also operate on battery power. This allows for measurements to be taken anywhere, even in places lacking a power source. The instrument can also be placed directly on the beaker, making it possible for quick and simple measurements to be taken. No need to establish a specific location for measurement — with VISCO, you can take measurements anytime, anywhere.

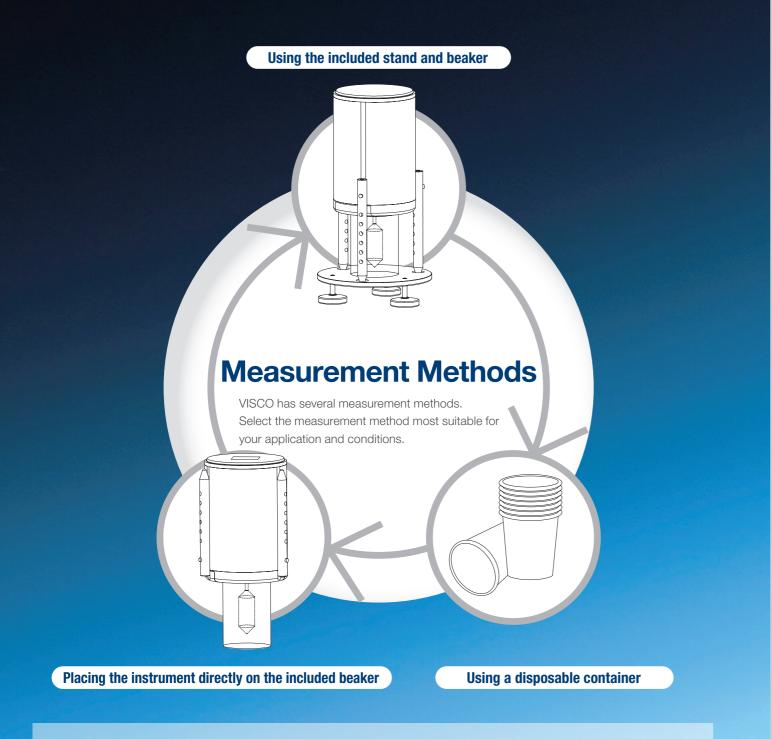

#### VISCO Package

#### Package A

Using disposable containers eliminates the hassle of cleaning after measurement. Package A, which includes a specialized adapter for use with disposable containers such as paper cups, is available at ATAGO.

• Cat.No.6810

 VISCO™ (main unit) · Cup Adapter (with 100pcs cups\*): RE-78141

- Cat.No.6830
- VISCO<sup>™</sup>-895 (main unit) · Cup Adapter (with 100pcs cups\*): RE-78141
- \* 50pcs of paper cups and 50pcs of plastic cups are included.

#### Package B

Measurement of low viscosity (1 to 2,000 mPa·s) is possible. A package that comes with Ultra Low Adapter (ULA) for measuring low viscosity sample and VISCO (main unit) is available.

#### Cat.No.6811

 VISCO™ (main unit) · Ultra Low Adapter (ULA): RE-77120

#### Cat.No.6831

Cup adapter setup example

Cup

adapter

· VISCO™-895 (main unit) · Ultra Low Adapter (ULA): RE-77120



<b>VISCO APPLI</b>	CATION
Food and Beverage	
Beverages (e.g. juice, etc.)	Viscosity is critical parameter in bev product reaches consumers.
Milk	In regards to milk, aside from whole (skim) milk. In general, nonfat milk h
Tomato Juice/Purée	Tomato juice or purée must always Viscosity management is indispensa
Sauce (Worcestershire sauce, thicker Worcester sauce, bork cutlet sauce, etc.)	There are many kinds of sauce. The thicker Worcester sauce and pork of determined by JAS.
Ketchup	Ketchup, a pseudoplastic fluid, is cl upside-down. Applying a bit of force for reacting differently at varying ten
Mayonnaise	Mayonnaise also remains in its bott viscosity. The greater the force appl
Olive Oil	There are many vegetable based oil when force is applied). Olive oil is a
Honey	Honey is a Newtonian fluid. Its visco change in viscosity.
Jam	Imagine spreading jam on a piece of making jam spreadable. Managing t
Yogurt	Numerous factors throughout the m pasteurization and pH managemen
Butter/Margarine	Butter is a Bingham plastic (a type of applied, but applying force past a c
Japanese Curry (curry roux)	Thickened curry (roux) is quite main changes it into a more paste-like co stage of being sealed into a retort p amount to extrude every time the se
Gelatin/Agar	Viscosity measurements can be use if the gelatin or agar completely soli preventing measurements from beir
Household Essentials	
Toothpaste	Toothpaste with a paste-like consist It is important for toothpaste to be a

flattening.

everage manufacturing, from the production phase all the way until the

ble milk (3.25%), there is reduced fat milk (2%), low-fat milk (1%) and nonfat has the lowest viscosity.

s flow through the production line under a constant, homogenized state. sable to this process.

nese include (in ascending order of viscosity): Worcestershire sauce, cutlet sauce. In Japan, there are approximate levels or grades for viscosity

characterized by its propensity to remain in its bottle even when turned ce (squeezing) to the bottle causes the ketchup to flow out. It is also known mperatures.

ttle, even when turned sideways or upside-down and maintains high blied, the easier it will flow out and the viscosity will decrease.

oils that are Newtonian fluids (a fluid that does not change viscosity even a Newtonian fluid.

cosity is not affected by force and speed. Only temperature can cause a

of toast. The jam easily glides across the toast. Viscosity is a crucial factor in the viscosity can be quite difficult, as jam contains solids.

manufacturing process, such as how much fat is left in the yogurt, ent affect the final product and texture (viscosity).

of non-Newtonian fluid). It can not flow unless some degree of force is certain degree cause it to become more malleable in proportion to the force.

nstream in Japan. Thickened curry is made by applying heat to flour, which consistency, resulting in an increase in viscosity. Even in the final processing pouch, the curry roux must maintain the same viscosity to allow the same same amount of force is applied.

sed to check and manage the gelling process of gelatin or agar. However, lidifies during viscosity measurement, a spindle-shaped gap will form, ing taken.

istency is a Bingham plastic. It will not flow out unless the tube is squeezed. be at optimal viscosity. After applying the appropriate amount onto a toothbrush, toothpaste at just the right viscosity will break cleanly from the tube and retain its shape without

#### Shampoo and Coditioner

Cosmetics



Conditioner must have a particularly high viscosity, as it coats every single stand of hair. The components used in shampoo and conditioner may not mix well, but adding viscosity ensures that they are evenly maintained.

Viscosity measurements and research in the rheological properties of cosmetics is conducted in order to give even slightly viscous cosmetics a smooth, light, easily spreadable quality when applied.

#### Industrial / Chemical

industrial / Onernical	
Adhesives	There are many types of adhesives for different purposes and applications. For example, structural adhesives include general adhesive for temporary bonding, gap-filling adhesive and adhesive used for coating purposes. <b>Gap-filling adhesives:</b> properties, such as bond strength, leveling (smoothness) and ease/difficulty of flow are assessed through viscosity measurements.
Water glass	Water glass is also commonly known as, "Sodium silicate." It is often used in soaps and adhesives. It is also used in a wide variety of fields, such as engineering, paper manufacturing and pharmaceuticals. Water glass has an extraordinarily high viscosity.
Resins/Polymers	Various kinds of resins and polymers have appeared in recent years. In addition to ascertaining their properties, viscosity assessment is also an absolute must.
Photopolymers (used in 3D printing)	With the spread of 3D printers, photopolymers have suddenly gained traction in the global market. When photopolymers are exposed to light and heat, their viscosity increases and they harden.
Paint / Pigment / Varnish / Insulators	<ul> <li>Paint (brush application): good brushability and drip-resistant. Tends to have a low viscosity under a high-shear rate and high structural viscous properties under a low-shear rate.</li> <li>Spray painting (coating): types of spray painting include air spray painting, airless spray painting and electrostatic coating. Most air spray paints have the same structural viscous properties as paint for brush applications. Airless spray painting and electrostatic coating is utilized in applications such as the final coating in automobile painting. As such, great emphasis is placed on the appearance and finish of the paint. Compared to paint for brush applications and air spray paint, most airless spray paints have a low viscosity, with similar properties to Newtonian fluids. The leveling of the paint after it's applied affects the overall appearance of the finished products.</li> <li>Electrodeposition (electrophoretic deposition): used for applying coatings to complex shapes/objects. Through electrodeposition, a film of coating is created on the surface of the target object. The object undergoes a baking or drying process, which makes the surface become smooth. This (leveling) is an extremely critical part of the process. It is necessary to use paint with viscosity sensitive to temperature.</li> </ul>
Glycerin	Glycerin is highly viscous and it's a great moisturizer. It is often added to cosmetics to increase their moisturizing properties. It is also used in various food products as a thickening agent and as a humectant in pharmaceuticals.
Hydraulic Fluids	The viscosity of hydraulic fluid tends to decrease as the temperature rises. If the viscosity decreases too much, it may lead to deterioration in the lubricating properties of the fluid and adversely affect pump efficiency. In contrast, viscosity tends to increase as the temperature decreases, this can increase the usual amount of force needed to operate the pump, and may accelerate wear and tear. This is why you must select a hydraulic fluid with the optimal viscosity for your application.
Lubricant	The viscosity of lubricant is a vital element, of which the life span is affected by the wear and tear of machinery. Depending on the viscosity, if too much heat is generated, the life span of a lubricant may be reduced. Not only does this increase costs, it also shortens the life span of the machinery. This is why managing the viscosity of lubricant is of the utmost importance.
Cutting Fluids	A lot of heat is generated during machining processes. Cutting fluid has a low viscosity, which makes it well-suited for effectively dissipating heat. Low viscosity is also necessary for facilitating circulation and eliminating impurities. However, to ensure precision and control when spraying or coating, the viscosity must not be too low. Managing the viscosity of cutting oil is essential.
Slurry/Ceramic Slurry (ceramic mold casting)	"Slurry" indicates a fluid which contains solid, suspended granules. Due to its properties, if it not mixed, it may precipitate. Viscosity management is critical for ensuring a smooth transition.

# **Thick (viscous) Food Measurement**

### VISCO-Thickness Meter

In recent years, Japan has faced an ever increasing, aging population. "Longevity" isn't the only point of discussion — how to extend good health throughout life expectancy and what defines QOL (Quality of Life) have also become hot topics.

Food is of course fuel for life. We essentially "eat to live," but food holds greater significance. Food can also bring us joy as a source of pleasure in our lives. Food consistency is an important component that impacts the quality of life for individuals that have difficulty in chewing and swallowing (dysphagia). VISCO can assist in managing food consistency, thereby improving the quality of life.

### The Importance of Thickening Food

A common misconception found among dysphagia diets and care facility diet plans is thickening food/drink (by grinding, pureeing, mashing, chopping, or mincing) makes it easier to swallow.


In reality, thickening food reduces the flow of the substance, and on the contrary, it becomes rather difficult to swallow. If this is the case, then just why is it important to thicken food?

Individuals with dysphagia (an affliction characterized by the failure in the automatic process of closing the respiratory tract to allow food to pass through the esophagus) need more time when swallowing to allow food to properly pass through the esophagus. Thickening food and beverages is effective in preventing food from accidentally passing through the respiratory tract.

## Enhancing the Way Food is Experienced

In the past, dysphagia diet assessment was based solely on "hardness," but recently, it has become common to assess dysphagia diets based on 3 criteria: hardness, cohesion (ease of clump formation) and adhesion (smoothness). There was also a time when the focus was on determining the best way of intaking food to absorb the most nutrition. Nevertheless, if you don't look forward to your meals or derive any enjoyment from eating in the first place, you may gradually stop eating much of anything. Flavor isn't the only aspect to consider; food texture (mouthfeel) is also regarded with great importance. Until now, from a number of standpoints such as cost, space (installation) and operation, large sized viscosity measurement devices and expensive texture meters were utterly ill-suited for personal use and for small-scale care facilities. VISCO is compact, lightweight and portable, making it ideal for hospitals, care facilities or household use. It can quickly and easily measure and assess the textual properties (thickness) of foods for dysphagia diets and care facility diet plans. Using VISCO, anyone can look forward to savoring and enjoying safe, nutritious food.

#### **Dysphagia Diet Pyramid**

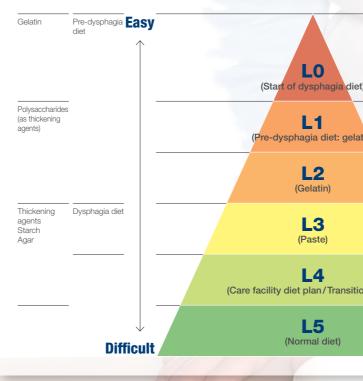

Omoyu gelatin (rice broth based gelatin), Kayu (rice gruel), Gelatin kayu (gelatin based rice gruel), Mochi (Japanese rice cake made of mochigome, a short-grain japonica glutinous rice), Negitoro (the fatty flesh of tuna minced to a paste and mixed with chopped green leaves of Welsh onions). Chawamushi (a cup-steamed egg custard hotchpotch), Foie gras (a food product made of the liver of a duck or goose that has been special) fatened), Mizuyoukan (a thick, soft jellied dessart made of red been paste), Koshian (boiled, mashed sweet red bean paste), Kabocha (Lapanese pumpkin; a variety of winter squash of the species (Cucurbita maxima), Shiitake (a variety of edible Japanese mushroom), Gomokumame (a Japanese dish consisting of cocked soybeans with various vegetables), Hijiki (a type of edible seaweed; Hizikia fusiforme)

Nutrition and Food	ł
BAD	GOOD
Intravenous (parenteral nutrition) Tube	e feeding (enteral feeding) Ingestion
Medical treatment for individuals unable to obtain adequate nutrients	Nutrients obtained through one's own accord (via ingestion).

	Dysphagia diet Main Foods			
	Grapes gelatin None desserts (Jell-O)		Uniform texture/ thickness	Start diet
	Omoyu gelatin	Negitoro, Chawanmushi		Dysphagia diet1
	Omoyu gelatin	Foie gras mousse		Dysphagia diet2
	Kayu, Gelatin kayu	Mizuyoukan, Egg dishes	Varying texture/ thickness	Dysphagia diet3
I diet)	Cooked rice, Kayu	Koshian, Kabocha, Stewed tender food		Transitional diet (care facility diet plan)
	Cooked rice, Mochi	Shiitake, Bread roll, Gomokumame, Hijiki		Normal diet